



COURAGE + KHAZAKA electronic GmbH, Mathias-Brüggen-Str. 91

50829 Köln, Germany

Tel.: +49 221-956499-0, Fax: +49 221-956499-1

info@courage-khazaka.de, www.courage-khazaka.de

Z0380/Z0384 Corneofix® F 20 (angular):

20 x 19.5 mm, dispenser box with 100 pcs. / roll of 2000 pcs.

Z0391/Z0381 Corneofix® D 100 (22 mm Ø):

disc, 22 mm Ø, dispenser box with 500 pcs. / roll of 5000 pcs.

Z0382/Z0837 Corneofix® D 102 (30 mm Ø):

disc, 30 mm Ø, roll of 5000 pcs.

Z0388/Z0389 Corneofix® D 104 (14 mm Ø)

disc, 14 mm Ø, roll of 5000 pcs.

The disc-shaped product comes with a plastic tweezer



Description & Intended use

The sticky tapes collect dry skin scales (desquamation) from the surface. The more and thicker the flakes on the tape, the dryer the skin. Sufficiently moisturized skin shows a thin and even film of corneocytes.

Take a Corneofix® off the roll by the supplied tweezer. Press the sticking part of the tape to the skin so that it sticks to the skin evenly. **Important:** Please make sure that your hands are washed and oil-free before touching the tape. Fingerprints on the foil may influence the measurement. Never touch the sticking side of the tape!

Take the Corneofix® off the skin and perform your analyses.

Technical Data

Corneofix® D 100/102/104: transparent tape for the desquamation measurement

Dimensions: 22, 30 or 14 mm Ø, thickness: approx. 0.1 mm


Tapes are completely harmless for the skin

A safety data sheet of the material can be obtained on request.


Optional accessory

Pressure-Aid-Instrument for all desquamation foils

Since constant contact pressure of the tape on the skin enhances reproducibility of measurements, the tool is equipped with an integrated spring. When you press the tape onto the skin against the pressure of this spring, a small pin will move upwards through the instrument shaft. This allows you to precisely control the pressure applied with your finger on top of the tool. Increase pressure until you start to feel the pin on your fingertip. Do not increase pressure further, just hold it for 10 seconds.

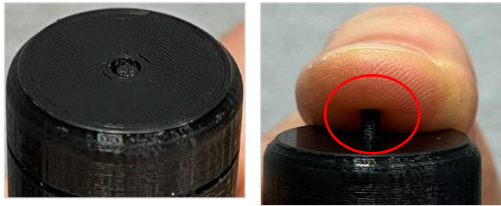

Correct pressure (2.2N) too much pressure applied

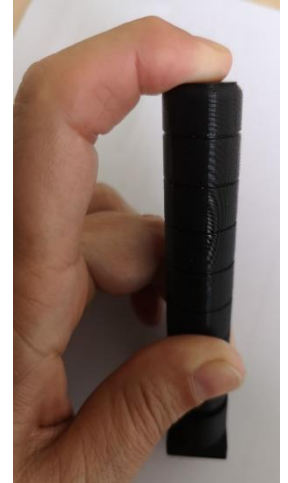

The Pressure-Aid-Instrument is not a measuring tool, but an aid for placing the Corneofix[®] tape evenly on the skin surface.

Technical Data

Pressure-Aid-Instrument:

Length: approx. 10 cm, weight: approx. 28 g, applied pressure: 2.2 N ~200 g/cm² (due to the different sizes of the pressure surface, the pressure per cm² can vary within a certain tolerance).

- Z0385** Pressure surface angular, 20 x 20 mm
- Z0385S** Pressure surface round, Ø 16 mm
- Z0385L** Pressure surface round, Ø 24 mm
- Z0385XL** Pressure surface round, Ø 32 mm



Z0385



Z0385S



Z0385L



Z0385XL

More questions? Please contact us.